

December 2024

These are the instructions for upgrading SAll software from a download link. PC-SAM software supports all of SA Instruments' single channel monitoring and gating systems. The current version of PC-SAM is 11.12. It has SnapView embedded. This version of PC-SAM supports numerous options including the Fluid Heater Systems, the Large Animal Capnograph and NIBP options. It also applies improvements controlling on animal temperature for the Fluid Heater Systems, to the Model 1030 internal battery charger and improvements in temperature measurement accuracy and fixes known bugs. Some of the windows are movable, others open away from displayed data.

PC-SAM.exe, TrendMap.exe and several support files reside in the directory "c:\ProgramFiles\PC-SAM".

To upgrade to the new versions, do the following:

1. Navigate to c:\ProgramFiles\PC-SAM or c:\ProgramFiles(x86)\PC-SAM
2. Copy and paste the PC-SAM folder in the ProgramFiles folder. This creates a backup of your previous version of PC-SAM.
3. Copy all files from the upgrade download into the PC-SAM folder. Answer yes to overwrite files.

Execute PC-SAM by clicking on the PC-SAM icon on the desktop. Enter the serial number from the bottom of your ERT Control/Gating Module, Control/Gating Module or Model 1025T. Navigate to PC-SAM's main display. PC-SAM revision 11.12 should be reported in the SYSINFO window.

If you get an error message when you try to execute one of the files that says "...you may not have authorization to execute this file...", do the following:

1. Navigate to the new .exe file,
2. Right click on the file name
3. Select Properties
4. Open the General Tab Page
5. At the bottom of the page, if it reads "This file came from another computer and might be blocked to help protect this computer" click the "unblock" button.

Problems contact SAll 631.689.9408.or service@i4sa.com.